
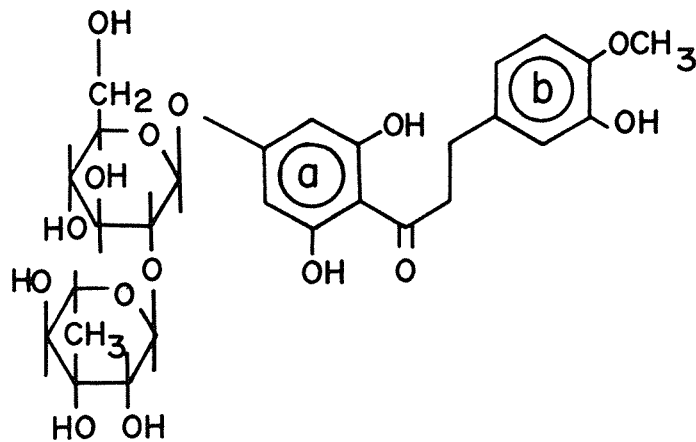


# CV NOTES

## NEOHESPERIDIN DIHYDROCHALCONE

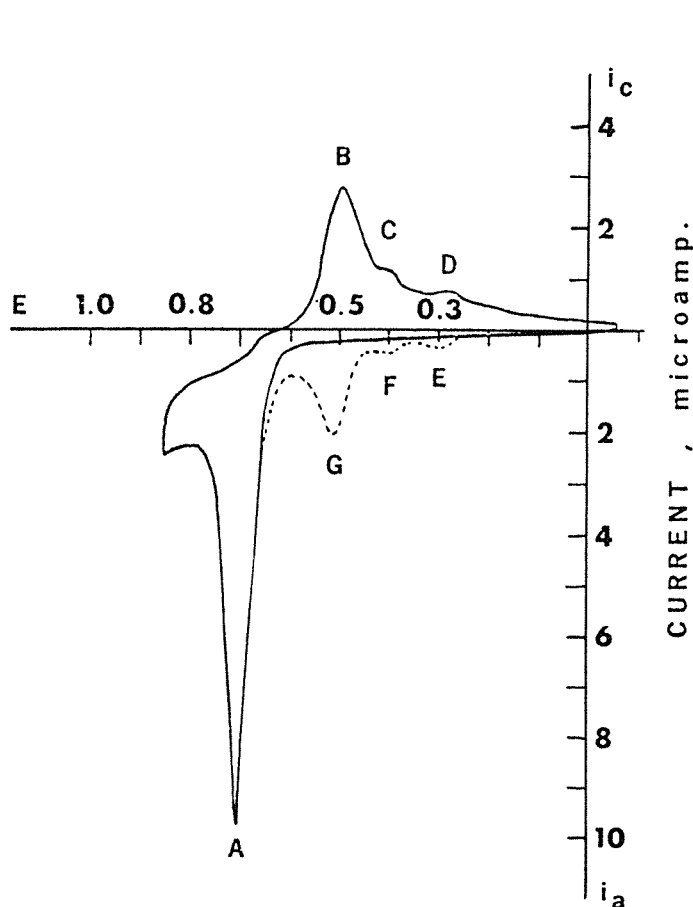

SAMPLE: Neohesperidin  
Dihydrochalcone

MEDIUM: 4 M HClO<sub>4</sub>

CONC: 2 mg/20 ml

RATE: 300 mV/s

ETRODE: CPE

REF: RE-1, Ag/AgCl

MODEL: CV-1A

A: Two electron oxidation of the B ring producing the o-quinone and methanol

B: Reduction of the B-ring quinone formed at A

C: unknown

D: unknown

E: unknown

F: unknown

G: Second scan oxidation of the B-ring catechol formed via wave B

2,3',6-trihydroxy-4-O-B-neohesperidosyl-4'-methoxy-dihydrochalcone is a natural product with a very sweet taste. It illustrates that electrochemistry can be useful for study of some rather complex phenolic natural products.

Submitted by Dr. Thomas Kenyhercz, Purdue University



2701 Kent Ave  
West Lafayette  
Indiana 47906