

3-Syringe Bracket Parts List (MD-1002)

<i>Part Number</i>	<i>Description</i>	<i>Quantity</i>
MF-3050	Bee 3-Syringe Clamp	1
MF-3051	Bee Carriage Extender Bar	1
MF-3052	Bee 3-Syringe Holder	1
MR-3348	4-40 x ¼ Flat Head Machine Screw	2

Instructions for Assembly of 3-Syringe Bracket

1. Refer to the figure below. Unscrew the thumbscrew from the 1-syringe holder on the Baby Bee pump. Remove the spring and the 1-syringe clamp from the thumbscrew. Insert the thumbscrew through the hole in the 3-syringe clamp (MF-3050) then through the spring. Screw the thumbscrew into the thread on the 3-syringe holder (MF-3052).
2. Remove the 1-syringe holder from the Baby Bee pump by unscrewing the four screws that attach it to the pump cover. Place the 3-syringe mount on the pump cover with the thumbscrew aligned over the corresponding hole in the cover. Attach the 3-syringe mount to the pump cover using the four screws from the 1-syringe mount.
3. Screw the carriage extender bar to the front of the carriage block using the two screws (MR-3348) supplied.

