

Dental Acrylic Cement

MD-1300

Instructions: Mix powder and liquid in a 3-4:1 ratio. Build the cement up around the groove in the guide cannula and the bone anchor screws. Apply with a disposable syringe, poly pipette or spatula.

Note: This product replaces the Sevriton dental acrylic cement previously sold as BASi part number MF-5181. Sevriton was discontinued by its manufacturer and is no longer available. BASi conducted an extensive replacement search and from the standpoint of cure time, exothermy, price and packaging, we think the enclosed cement is an appropriate replacement.

Please address comments to:

Microdialysis Group	765-463-4527
Bioanalytical Systems, Inc.	765-497-1101 fax
2701 Kent Avenue	
West Lafayette, IN 47906 USA	